

PRODUCTS FOR GROWTH PROMOTION TESTING

The broadest collection of QC microorganism products for Growth Promotion Testing

Microbiologics



Labcare
de Colombia



	EZ-Accu Shot™ Select	EZ-Accu Shot™	EZ-CFU™ One Step	EZ-CFU™
Available Strains	5 compendial strains plus <i>E. coli</i> in one convenient kit	40+ licensed, traceable strains	20+ licensed, traceable strains	20+ licensed, traceable strains
Number of Tests	10 tests per vial 10 tests for each organism per kit	10 tests per vial 50 tests per kit	19 tests per vial 190 tests per kit	90+ tests per vial 900+ tests per kit
Hydration Fluid (Included in Kit)	Equilibrate to room temp	Equilibrate to room temp	Pre-warm Hydrating fluid 30 minutes	Pre-warm Hydrating fluid 30 minutes
Lyophilized Microorganism Pellets	Use 1 pellet per suspension to obtain 10-100 CFU / 0.1 ml	Use 1 pellet per suspension to obtain 10-100 CFU / 0.1 ml	Use 2 pellets per suspension to obtain 10-100 CFU / 0.1 ml	Use 2 pellets per suspension to obtain 10-100 CFU / 0.1 ml
Resuscitation of Microorganisms	Quick-dissolve pellet, no pre-incubation	Quick-dissolve pellet, no pre-incubation	Incubate suspension for 30 minutes before inoculation	Incubate suspension for 30 minutes before inoculation
Dilutions	No dilution step	No dilution step	No dilution step	One log dilution required
Stability of Hydrated Suspension	8 hour stability	8 hour stability*	8 hour stability*	30 minute stability

* Please see instructions for use for 8 hours stability exceptions

Choose from 4 unique products that meet your specific needs – all designed to deliver 10-100 CFU as required by the USP/Ph. Eur./JP.